

## **Fluid and Filters**

### **Cooling Fluid**

Check the cooling fluid level daily, top up as necessary. Weekly, empty the cooling fluid bottle and refill with clean fluid.

### **Fluid Filter**

Check the fluid filter every 2 months. This is done by simply lifting the suction pipe out of the reservoir bottle, unscrewing the original filter and fitting a new one then replace the suction pipe in the reservoir bottle.

### **Air Filter**

Examine and clean the air filter regularly. Do not operate the instrument without the air filter in place, foreign bodies can be drawn into the cooling system and impair cooling efficiency. It is good practice to have a spare clean filter to replace the dirty one. The element should be washed and thoroughly dried.