

Loading Printer Paper

1/ Press down the printer cover release button located immediately in front of the printer cover, then lift up on the printer cover to open. Release the printer cover release button.


2/ Unroll about 15 cm of paper from the paper roll.

3/ Remove the old roll and insert the roller into the new paper roll.

4/ Place the paper roll and roller into the V-shaped roller supports. Make sure the paper roll is feeding from the bottom of the roll. When both ends of the roller are in proper position, the paper roll should be level. If the paper roll is not level, carefully reseal the roller between the roller guides.

5/ Hold the paper down onto the surface of the Cryoscope and centre the paper on the printer tear bar. When the paper is centred, gently close the printer cover. Press on the top of the printer cover to make sure it is firmly latched with the printer mechanism.

6/ The printer paper is loaded and ready for operation.

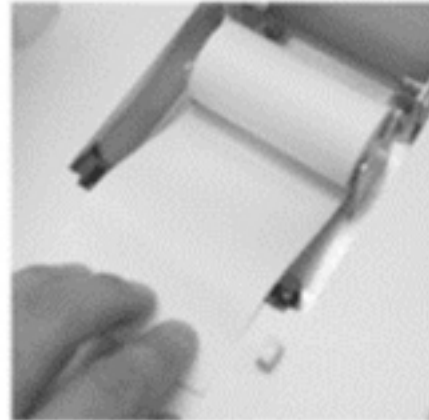

Caution

Do not attempt to pull the paper through the printer, as this may damage the mechanism. Either release the printer cover by fully depressing the printer cover release button and lifting up on the cover, or use the Feed button on the keypad to advance the paper.

