



Absorbance difference between Colour and Bitterness cuvettes for the BeerLab

1. From the settings menu, select 'Reading Test'

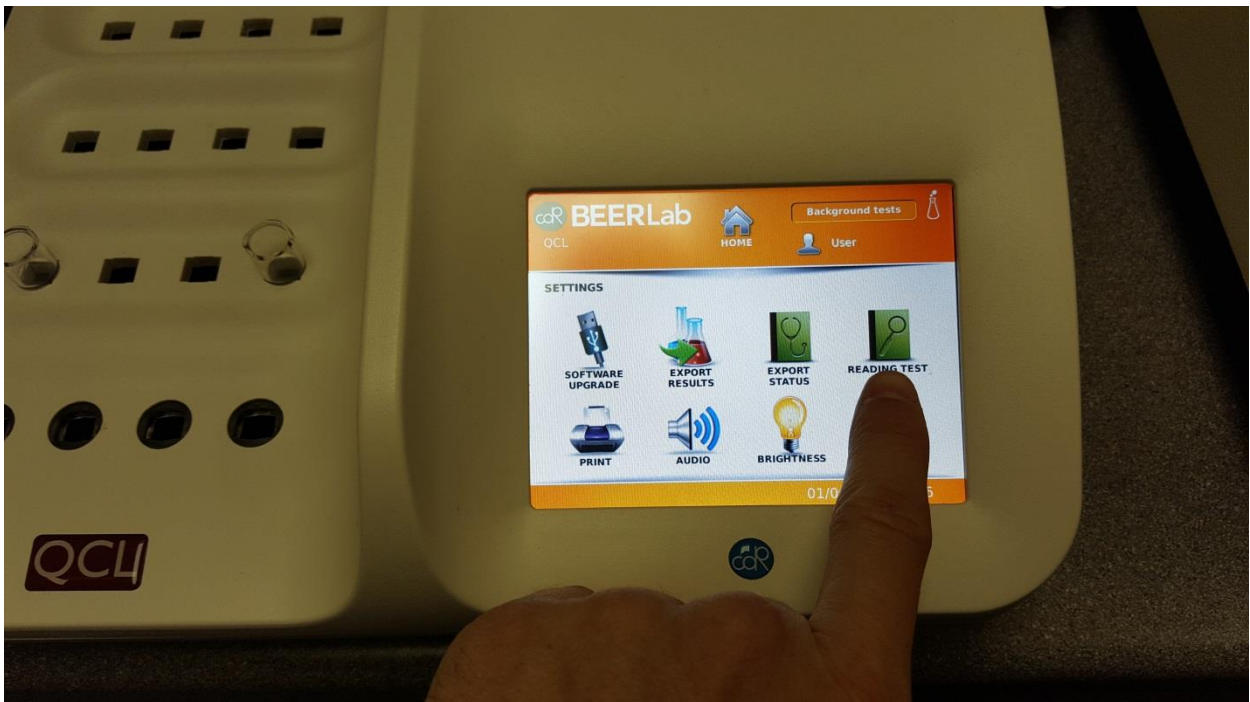

2. On the 'Reading Test' screen select cuvette LED I as shown below

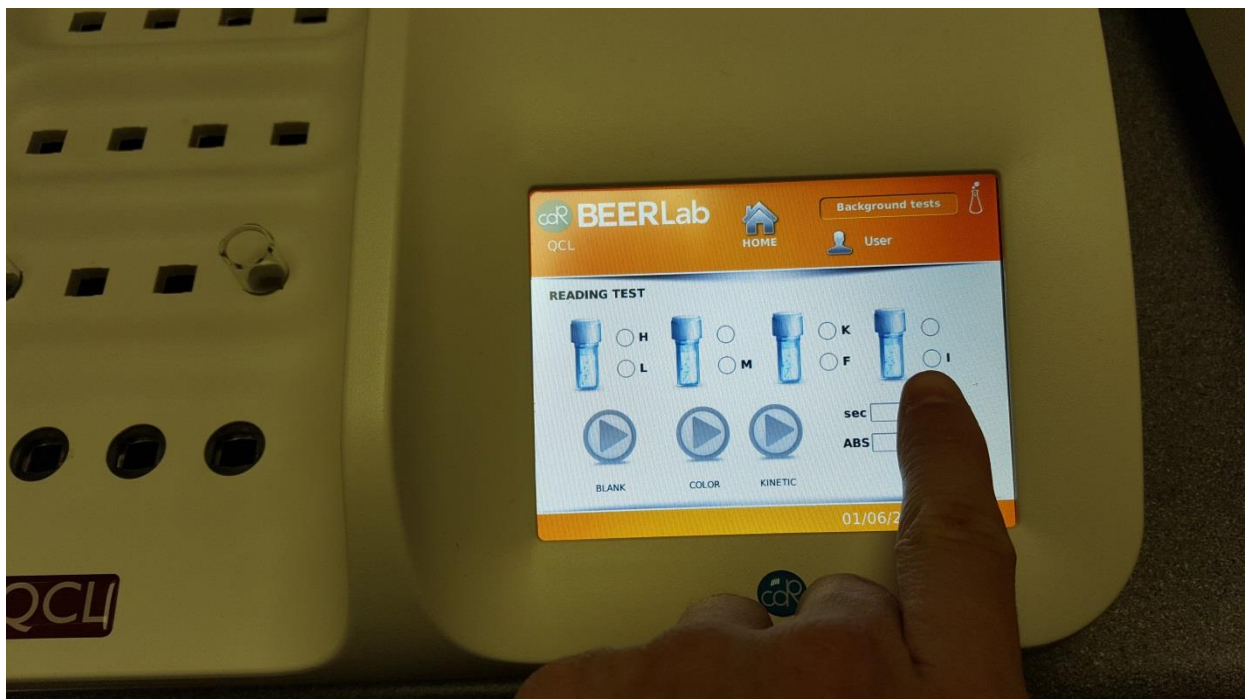

3. Press 'Blank' with nothing in the reading cell



4. Place a cuvette in the reading cell marked with blue light and press 'colour' – The cuvette below is a bitterness cuvette.



5. The reading test was repeated with a colour cuvette, showing a near 10-fold increase in absorbance



In summary, bitterness cuvettes have an absorbance around 1.0000 to 1.5000 and colour cuvettes have an absorbance around 11.0000.

If your colour and bitterness cuvettes have been mixed up, this is a quick and easy method of distinguishing the two.